



PeliKine Compactä human IL-8 ELISA kit

288 tests

An enzyme immunoassay for the quantitative determination
of human Interleukin 8

PRODUCT INFORMATION

Sanquin Reagents

Plesmanlaan 125
1066 CX Amsterdam
The Netherlands

Cat.No. M1918

For The Netherlands:

Tel. 020.512 3246

Fax. 020.512 3570

For other countries:

contact your local distributor.

Protocol summary and checklist PeliKine compactă human IL-8 ELISA kit

Day 0:

- Bring coating antibody to room temperature (18-25 °C).
- Prepare coating buffer.
- Dilute coating antibody 1:100 in coating buffer, add 100 µl to all wells, cover the plate(s) and incubate overnight at room temperature.

Day 1:

- Bring all reagents, with the exception of streptavidin-HRP, to room temperature.
- Prepare blocking buffer
- Wash the plate(s) five times with PBS.
- Add 200 µl blocking buffer to all wells and incubate for one hour at room temperature.
- Prepare dilution buffer, standard and sample dilutions.
- Prepare washing buffer.
- Wash the plate(s) five times with washing buffer.
- Leaving the substrate blank wells empty, add 100 µl of standard and sample dilutions to the appropriate wells, cover the plate(s) and incubate for one hour at room temperature.
- Dilute biotinylated antibody 1:100 in dilution buffer.
- Wash the plate(s) five times with washing buffer.
- Leaving the substrate blank wells empty, add 100 µl of the diluted biotinylated antibody to all wells, cover the plate(s) and incubate for one hour at room temperature.
- Dilute the streptavidin-HRP conjugate 1:10,000 in dilution buffer.
- Wash the plate(s) five times with washing buffer.
- Leaving the substrate blank wells empty, add 100 µl of the streptavidin-HRP conjugate to all wells, cover plate(s) and incubate for 30 minutes at room temperature.
- Just before use, prepare substrate solution.
- Wash the plate(s) five times with washing buffer.
- Add 100 µl substrate solution to all wells, including the substrate blank wells, and incubate for 30 minutes at room temperature in the dark.
- Add 100 µl stop solution to all wells and read the plate at 450 nm.
- Calculate the amount of IL-8 in the samples.

	1	2	3	4	5	6	7	8	9	10	11	12
A	S1	S1										
B	S2	S2										
C	S3	S3										
D	S4	S4										
E	S5	S5										
F	S6	S6										
G	S7	S7										
H	S8	S8								B	B	

Plate plan proposed for the PeliKine compact™ human IL-8 ELISA kit:

Key: B: substrate blank S1-S8: IL-8 standards 0 - 240 pg/ml Empty: samples

I. INTRODUCTION

Interleukin 8 (IL-8) is a cytokine with multifunctional actions. Several substances originally described for their biological activities have been identified as IL-8; monocyte-derived neutrophil chemotactic factor (MDNCF), neutrophil-activating peptide (NAP), neutrophil-activating factor (NAF), granulocyte chemotactic protein (GCP), T-lymphocyte chemotactic factor (TCF), leukocyte adhesion inhibitor (LAI). IL-8 was isolated first from stimulated leukocytes, but the molecule can be produced by a wide variety of cell types in response to cytokine inducers. In view of the increasing number of newly discovered chemotactic proteins, that are structurally related to IL-8, these cytokines are now designated 'chemokines'. The hallmark for this family of proinflammatory proteins is the conservation of four cysteine residues that are important for the tertiary structure. The chemokines can be divided into two subfamilies depending on whether the first two cysteines are adjacent (C-C chemokines) or not (C-X-C chemokines). IL-8 is a member of the second subfamily.

Elevated levels of IL-8 have been after sublethal endotoxaemia, septic shock, microbial infection of the amniotic cavity, Jarish-Herxheimer reaction of relapsing fever, infectious diseases of the central nervous system, acute pancreatitis, ulcerative colitis, empyaema, haemolytic uraemic syndrome, meningococcal disease, gastric infection, pertussis, and peritonitis.

Bioassays for the quantification of IL-8, based on the ability to chemoattract neutrophils in vitro, have been used for several years. These assays, although sensitive, are time consuming and susceptible to interference by other substances.

This PeliKine compact™ human IL-8 kit has been developed for faster, more reproducible and specific quantification of human IL-8 (huIL-8) in plasma and other body fluids, as well as in cell-culture supernatant.

II. PRINCIPLE OF THE TEST

The PeliKine compact™ human IL-8 ELISA kit is a "sandwich-type" of enzyme immunoassay in which a monoclonal anti-huIL-8 antibody is bound onto polystyrene microtiter wells. Human IL-8, present in a measured volume of sample or standard is captured by the antibody on the microtiter plate, and non-bound material is removed by washing. Subsequently, a biotinylated sheep antibody to huIL-8 is added. This antibody binds to the huIL-8-antibody complex present in the microtiter well. Excess biotinylated antibody is removed by washing, followed by addition of horseradish peroxidase (HRP) conjugated streptavidin, which binds onto the biotinylated side of the huIL-8 sandwich. After removal of non-bound HRP conjugate by washing, a substrate solution is added to the wells. A coloured product is formed in proportion to the amount of huIL-8 present in the sample or standard. After the reaction has been terminated by the addition of a stop solution, absorbance is measured in a microtiter plate reader. From the absorbance of samples and those of a standard curve, the concentration of huIL-8 can be determined by interpolation with the standard curve.

III. STORAGE AND STABILITY

The PeliKine compact™ human IL-8 kit should be stored at -18°C to -32°C . The performance of the PeliKine compact™ human IL-8 kit is guaranteed until the expiration date shown on the case label.

IV. CONTENTS OF THE KIT


The PeliKine compact™ human IL-8 kit contains material sufficient for 288 tests, including standard curve samples. The reagents provided are:

Quantity	Kit component	Volume	Cap colour
1 vial	coating antibody 100-fold concentrated	375 µl	red
1 vial	blocking reagent 50-fold concentrated	2 ml	transparent
1 vial	IL-8 standard 10 ng/ml	200 µl	black
1 vial	biotinylated IL-8 antibody 100-fold concentrated	375 µl	yellow
1 vial	streptavidin-HRP conjugate 10,000-fold concentrated	20 µl	brown
1 bottle	dilution buffer 5-fold concentrated	60 ml	-
3 pcs	microtiter plates + lid	-	-
10 pcs	plate seals	-	-

V. PRECAUTIONS FOR USE

- 1) The PeliKine compact™ human IL-8 kit is intended *for research purposes only*.
- 2) Only use the reagents and microtiter plates supplied with the kit, do not mix reagents from different kit lots.
- 3) Handle all plasma and serum samples with care to prevent transmission of blood-borne infections.
- 4) Sodium azide inactivates HRP, do not use sodium azide-containing solutions, nor add sodium azide to the supplied materials.
- 5) All reagents contain thiomersal (0.001 % w/v) and may be toxic upon ingestion, inhalation or skin contact. Avoid contact of skin, eyes or clothing with dilution, washing or substrate buffer. In case of contact, wash skin or eyes with water and consult a physician.
- 6) The IL-8 standard contains human serum which has been tested for the markers of specific disease-transmitting agents in accordance with current EU guidelines to GMP and found to be nonreactive. Nevertheless the standard should be handled as potentially hazardous and capable of transmitting diseases.
- 7) Centrifuge all vials before use (1 minute 3000 x g)

VI. ADDITIONAL BUFFERS & SOLUTIONS REQUIRED

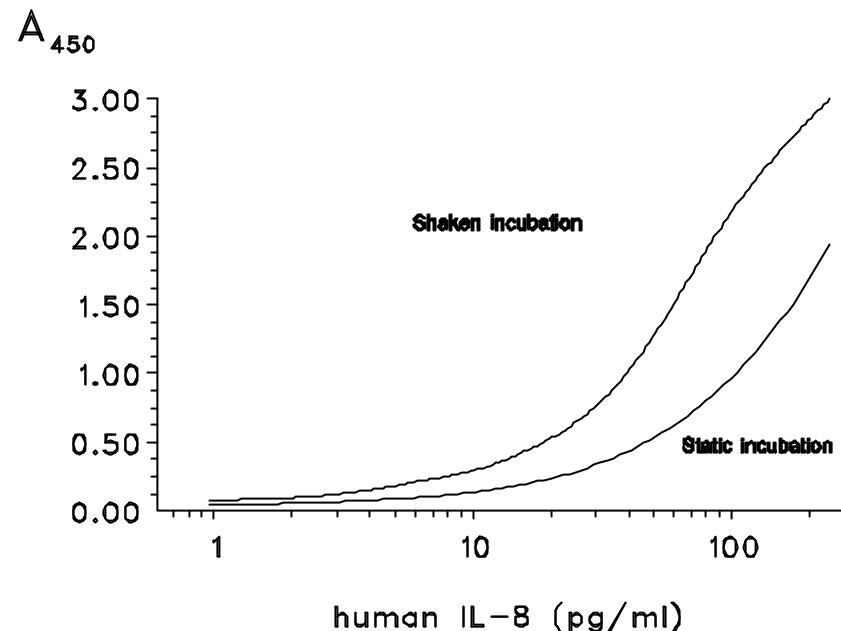

Typical standard curve for PeliKine compact™ human IL-8 ELISA kit
The assay is completed static or shaken at room temperature.

		STATIC INCUBATION	SHAKEN INCUBATION
		Calculated mean absorbance at 450 nm	
substrate blank		0	0
0	pg/ml	0.038	0.041
1	pg/ml	0.041	0.077
2.5	pg/ml	0.060	0.107
6.1	pg/ml	0.097	0.215
15.4	pg/ml	0.189	0.424
38.4	pg/ml	0.418	0.974
96	pg/ml	0.940	2.130
240	pg/ml	1.943	3.000

DO NOT USE THESE DATA TO CONSTRUCT A STANDARD CURVE FOR SAMPLE VALUE CALCULATIONS

14. STOP ENZYMATIC REACTION

Add 100 µl of stop solution to all wells.

After stopping the colour is stable for maximally 30 minutes.

15. PLATE READ-OUT

Read at 450 nm in an ELISA reader.

IX. RESULTS**Substrate blank**

- Record the absorbance at 450 nm for the substrate blank wells and average the duplicate values.

Standard curve

- Record the absorbance at 450 nm for each well containing standard and average the duplicate values.
- Calculate the net average absorbances by subtracting the average of the substrate blank wells.
- Plot the net average absorbances (ordinate) versus the IL-8 concentration in pg/ml (abscissa) on log-linear paper and draw the best fitting curve. An example of a standard curve is given on the next page.

Samples

- Record the absorbance at 450 nm for each standard well, and average the duplicate values.
- Calculate the net average absorbances by subtracting the average of the substrate blank wells.
- Locate the net average absorbance value found for each sample on the vertical axis and follow a horizontal line intersecting the standard curve. At the point of intersection, read the IL-8 concentration (pg/ml) from the horizontal axis. Multiply the obtained IL-8 concentration with the dilution factor of the sample and record this figure.

X. INCREASED SENSITIVITY

The assay sensitivity can be increased by a small adaptation of the incubation methodology. Just follow all the instructions as stated in the assay procedure (chapter VIII), but incubate at room temperature (18-25°C) on a horizontal plate shaker at 700 ± 100 rpm. All incubations, including the enzymatic colour development, have to be completed on the shaker, in the same time as stated in the static assay procedure. This will result in an increase in assay sensitivity, with small effects on the background levels (see figure next page).

Coating buffer: 0.1 M Carbonate/bicarbonate buffer pH 9.6

Solution A: 1.24 g $\text{Na}_2\text{CO}_3 \cdot \text{H}_2\text{O}$ in 100 ml distilled water

Solution B: 1.68 g NaHCO_3 in 200 ml distilled water

Take 70 ml of solution A, and add solution B until the pH is 9.6 (approximately 175 ml of solution B required)

The prepared buffer can be stored up to one week at 2-8°C.

PBS stock solution [20 x]: 0.2 M Phosphate Buffered Saline (PBS)

Dissolve	32 g	$\text{Na}_2\text{HPO}_4 \cdot 2\text{H}_2\text{O}$
	6 g	$\text{NaH}_2\text{PO}_4 \cdot 2\text{H}_2\text{O}$
	164 g	NaCl

in 900 ml distilled water

(intensive stirring and some heating will speed dissolution).

Bring the temperature of the solution back to room temperature (18-25 °C) and check pH; if necessary adjust pH to 6.8 - 6.9 with concentrated HCl or NaOH, and add distilled water to a volume of 1 liter (when diluted 20 times the obtained buffer will have a pH of 7.2 - 7.4).

Add 20 mg thiomersal as preservative. Do not use sodium azide (NaN_3) since this preservative reduces the quality of the enzymatic label.

The prepared buffer can be stored up to three months at 2-8 °C.

Note: in the concentrated buffer salt crystals may appear when stored at 28°C. Before preparing the working-strength buffer, first warm the concentrated buffer BRIEFLY to 37°C to dissolve the precipitate.

Washing buffer: PBS with 0.005 % TWEEN 20

Make 1 liter of working-strength PBS by diluting the PBS stock solution (see above) 20-fold with distilled water.

Add 50 µl TWEEN 20.

The prepared buffer can be stored up to one month at 2-8 °C.

Substrate buffer: 0.11 M acetate buffer pH 5.5

Dissolve 15.0 g sodium-acetate ($\text{CH}_3\text{COONa} \cdot 3\text{H}_2\text{O}$) in 800 ml distilled water.

Adjust pH to 5.5 with glacial acetic acid, add distilled water to a volume of 1 liter.

Do not add any preservative (e.g. merthiolate, sodium azide) since this may affect the quality of the enzymatic colour development.

The prepared buffer can be stored up to two weeks at 2-8 °C.

3,5,3',5'-tetramethylbenzidine (TMB) stock solution: 6 mg/ml TMB in DMSO

Dissolve 30 mg 3,5,3',5'-tetramethylbenzidine (TMB) in 5 ml dimethylsulfoxide (DMSO).

The prepared stock solution can be stored up to 1 month **at room temperature (18-25°C)** and **protected against light**.

Hydrogen peroxide stock solution: 3% H_2O_2 solution in distilled water.

The prepared stock solution can be stored up to one month at 2-8 °C.

Substrate solution

For each plate mix the following reagents:

- 12 ml substrate buffer
- 200 μl TMB stock solution
- 12 μl H_2O_2 stock solution

The substrate solution should be prepared just before use and has to be at room temperature (18-25°C) for optimal reproducible results.

Stop solution: 1.8 M H_2SO_4 solution in distilled water.**10. THIRD WASH STEP**

Aspirate supernatant from wells and wash the microtiter plate(s) as described in point 6 above.

11. THIRD INCUBATION STEP**Streptavidin-HRP conjugate**

The kit contains one brown capped vial of concentrated streptavidin-HRP conjugate, which must be stored at -18°C to -32°C to maintain maximal stability. The contents of the vial will not be frozen at this temperature.

Per microtiter plate, add 3 μl streptavidin-HRP conjugate to 30 ml of working-strength dilution buffer just before use. **Do not prepare in advance of assay.**

Leaving the substrate blank wells empty, add 100 μl of streptavidin-HRP conjugate to all wells.

Cover the microtiter plate(s) with adhesive seal, agitate the microtiter plate by tapping the edge of the frame for a few seconds to mix contents of each well and **incubate for 30 minutes at room temperature (18-25°C)**.

12. FOURTH WASH STEP

Aspirate supernatant from wells and wash the microtiter plate(s) as described in point 6 above.

13. FOURTH INCUBATION STEP**Enzymatic colour development**

Approximately 10 minutes before use, prepare the substrate solution as described on page 4 of this leaflet.

The substrate solution should be at room temperature (18-25°C) for optimal reproducible results.

Add 100 μl of substrate solution to all wells, **including the substrate blank wells**.

Cover microtiter plate(s) with lid, gently agitate the microtiter plate by tapping the edge of the frame for a few seconds to mix contents of each well and **incubate for 30 minutes at room temperature (18-25 °C) in the dark**.

do not cover the plate with aluminium foil.

Note: The speed of enzymatic colour development is influenced by many factors including temperature and quality of the used TMB.

6. FIRST WASH STEP

Prepare washing buffer as described on page 3 of this leaflet.

Wash the required microtiter plates five times with washing buffer in a plate washer.
In case of manual washing, completely fill the wells (> 300 µl) with washing buffer and aspirate, repeat this four times. After the final aspiration the wells should be dry.

7. FIRST INCUBATION STEP**Standards and samples**

Leaving the substrate blank wells empty, transfer 100 µl of the prepared standards and samples in duplicate into the appropriate wells (see recommended plate plan).
Cover plate(s) with adhesive seal, gently agitate the microtiter plate by tapping the edge of the plate for a few seconds to mix contents of each well and **incubate for 1 hour at room temperature (18-25°C)**.

8. SECOND WASH STEP

Aspirate supernatant from wells and wash the microtiter plate(s) as described in point 6 above.

9. SECOND INCUBATION STEP**biotinylated antibody**

The kit contains one yellow-capped vial with concentrated antibody-biotin conjugate.

Per microtiter plate, add 120 µl biotinylated antibody to 12 ml working-strength dilution buffer just before use.

Leaving the substrate blank wells empty, add 100 µl of diluted biotinylated antibody to all wells.

Cover plate(s) with adhesive seal, gently agitate the microtiter plate by tapping the edge of the plate for a few seconds to mix contents of each well and **incubate for 1 hour at room temperature (18-25°C)**.

VII. ADDITIONAL INFORMATION**Additional materials required**

- Pipetting devices for accurate delivery of 1-10 µl, 50 µl, 100 µl and 1 ml volumes.
- Beakers, flasks, cylinders necessary for preparation of reagents.
- Device for delivery of washing buffer (wash bottle / automated plate washer).
- Microtiter plate reader.

Sensitivity

MEAN calculated zero signal + 3 SD : 1 - 3 pg/ml (shake - static incubation)
2 x (MEAN calculated zero signal) : 4 - 8 pg/ml (shake - static incubation)

Note: the sensitivity is dependent of the type and quality of enzymatic substrate.

Expected values

IL-8 values in fresh serum and plasma samples of healthy individuals are below 10 pg/ml.

Specificity

No crossreactivity was observed with the following recombinant human proteins: IL-1 α, IL-1β, IL-2, IL-3, IL-4, IL-5, IL-6, IL-7, IL-9, IL-10, IL-11, Macrophage Colony Stimulating Factor (M-CSF), Granulocyte Colony Stimulating Factor (G-CSF), Granulocyte/Macrophage Colony Stimulating Factor (GM-CSF), Leukaemia Inhibitory Factor (LIF), RANTES, Stem Cell Factor/ Mast Cell Factor (SCF/MCF), Transforming Growth Factor β-1 (TGFβ-1), Tumour Necrosis Factor α (TNFα), Tumour Necrosis Factor β (TNFβ/Lymphotoxin), and Interferon γ (IFNγ).

REFERENCES

- Aggerwal, B. *et al* (1992) Human Cytokines ISBN 0-86542-183-8
 Baggiolini, M. *et al* (1993) Immunol.Today **14**: 24
 Broaddus *et al* (1992) Am.Rev.Respir.Dis. **146**: 825
 Crabtree *et al* (1993) Scand.J.Immunol. **37**: 65
 Friedland J.S. *et al* (1992) Infect.Immun. **60**: 2402
 Gross *et al* (1992) Eur.J.Clin.Invest. **22**: 200
 Hack, C.E. *et al* (1992) Infect.Immun. **60**: 25
 Halstensen *et al* (1993) J.Infect.Dis. **167**: 471
 Lin *et al* (1993) Nephron **63**: 404
 Martich *et al* (1991) J.Exp.Med. **173**: 1021
 Matsushima, K. *et al* (1989) Cytokine **1**: 2
 Negussie *et al* (1992) J.Exp.Med. **175**: 1207
 Sheron and Williams (1992) Clin.Exp.Immunol. **89**: 100
 Thomson, A.W. (1994) The cytokine handbook. ISBN 0-12-689661-5
 Torre *et al* (1992) Am.J.Dis.Child. **147**: 27
 Van Deventer *et al* (1993) J.Inf.Dis. **167**: 461
 Van Meir *et al* (1992) Cancer Res. **52**: 4297
 Van Zee *et al* (1991) J.Immunol. **146**: 3478

VIII. ASSAY PROCEDURE

1. **BRING ALL REAGENTS TO ROOM TEMPERATURE (18-25°C)**, with the exception of the streptavidin-HRP conjugate which has to be kept at -18 °C to -32 °C to ensure stability. Centrifuge all vials before use (1 minute 3000 x g).

For your convenience an easy-reference manual with check list and plate plan are available on the last pages of this leaflet.

2. DILUTION BUFFER

The kit contains one bottle with 5-fold concentrated dilution buffer.

Calculate the quantity of dilution buffer required (approximately 15 ml undiluted buffer per microtiter plate) and prepare a working-strength solution by diluting the opalescent concentrated buffer 5 times in distilled water before use. The working-strength dilution buffer can be stored for up to one week at 2-8 °C.

For optimal assay results, dilute samples and standard in working-strength dilution buffer.

3. MICROTITER PLATES

Coating antibody

Coating

The kit contains three microtiter plates for 96 tests each, including the standard curve samples.

Prepare coating buffer as described on page 3 of the information leaflet.

Per microtiter plate, add 120 µl of coating antibody (red-capped vial) to 12 ml coating buffer.

Add 100 µl to all wells, cover microtiter plate(s) with lid and **incubate overnight at room temperature (18-25 °C)**.

Washing procedure

Prepare working-strength PBS (1:20 dilution of stock PBS as described on page 3 of the information leaflet).

Aspirate supernatants from wells and completely fill the wells (> 300 µl) with working-strength PBS and aspirate. Repeat this four times, after the final aspiration the wells should be dry.

Blocking procedure

The kit contains one transparent-capped vial with 2 ml blocking reagent.

Prepare blocking buffer by adding 500 µl blocking reagent to 25 ml working-strength PBS (1:20 dilution of stock PBS as described on page 3 of the information leaflet).

Add 200 µl blocking buffer to all wells, cover microtiter plate(s) with adhesive seal, gently agitate the microtiter plate by tapping the edge of the plate for a few seconds to mix contents of each well and **incubate for 1 hour at room temperature (18-25 °C)**.

4. IL-8 STANDARD

Standard curve preparation

A natural huIL-8 standard has been calibrated against the WHO Interim International Standard (IL-8 89/520; National Institute for Biological Standards and Control, Potters Bar, Hertfordshire, U.K. 1 WHO Unit = 10 ng IL-8). In former CLB IL-8 reagents sets [Batch IL8-CK0001 to IL8-CK0004 and 1918-00-05] 1 pg IL-8 standard is comparable with 0.56 pg of the WHO standard).

The kit contains one black-capped vial with 10 ng/ml natural IL-8.

Label 7 tubes, one tube for each dilutions: 240, 96, 38.4, 15.4, 6.1, 2.5 and 1 pg/ml.

Pipette 610 µl of working-strength dilution buffer into the tube labelled 240 pg/ml and 300 µl of working-strength dilution buffer into the other tubes.

Transfer 15 µl of the IL-8 standard (10 ng/ml) into the first tube labelled 240 pg/ml, mix well and transfer 200 µl of this dilution into the second tube labelled 96 pg/ml.

Repeat the serial dilutions six more times by adding 200 µl of the previous tube of diluted standard to the 300 µl of dilution buffer.

The standard curve will contain 240, 96, 38.4, 15.4, 6.1, 2.5, 1 and 0 pg/ml (dilution buffer).

It is recommended to prepare two separate series for each assay.

Avoid repeated freeze-thawing of the standard, although experimental data have shown that up to 3 freeze-thaw cycles have no effect on the IL-8 levels of the standard.

5. SAMPLES

Serum, EDTA-anti-coagulated plasmas, and culture fluids are suitable for use in the assay (**caution:** separate plasma/serum and blood cells within 4 hours after collection, non-separated samples must be kept on temperatures from 2 to 8°C). Do not use grossly haemolyzed or lipemic specimens. If samples are to be run within 24 hours, they may be stored at 2-8°C; otherwise samples should be stored frozen (<-18°C).

Up to 3 freeze-thaw cycles have no effect on the IL-8 levels of serum or plasma samples. Nonetheless, excessive freeze-thaw cycles should be avoided. Prior to the assay, frozen samples should be thawed **as quickly as possible** in a 37°C waterbath and then brought to room temperature (18-25°C).

It is recommended to dilute the test samples at least 1:2 in working-strength dilution buffer. If high levels of IL-8 (>300 pg/ml) are expected in the test samples, additional dilutions of sample i.e. 1:10 and 1:50 should also be prepared.

